



**BIOL@GLX<sup>®</sup>**

**Biologix Group Limited**

**Address: No.2766 Ying Xiu Road High-Tech Industrial  
Development Zone Jinan, Shandong China**

**Phone: +86-531-67802668**

**Email: [Inquiry@BiologixGroup.com](mailto:Inquiry@BiologixGroup.com)**

**CKS-4008/CKS-5008**

**Multi-channel motorize decapper**

**Operation Manual**



**Biologix Group Limited**

## Foreword

Thank you for purchasing our Products: Multi-channel motorize decapper. This Manual for users contains function and operation of the Instrument. In order to use the instrument properly, please read this manual carefully before using the Instrument.

Multi-channel motorize decapper is used in the laboratory for all kinds of cryovial decapper or decapper . It is recognized as an efficient and excellent instrument in medical biochemistry analysis fields.

## Opening Check

Please check the Instrument and accessories with the packing list when you open the instrument packing case first time. If you find there is something wrong with the Instrument and the accessories ,please contact the vendor or the producer.

## Structure Description



### Multi-channel motorize decapper Operationinstructions:

A. Hold the decapper according to the picture and press the capper button and will make the capper action,align with the tube cap and press the decapper button and will make the decapper action.

B. Hold the decapper according to the picture and press the ejected button, The cap ejected will automatically .

C. If the decapper show the red indicator when operating , it indicates power is low. Please charge it in time. In order to better guarantee battery life and It must be recharged within 10 months,If you don't use the device for a long time.



### Performance parameter :

Capper/ Decapper speed	sec	4sec
Total weight	g	380g
Charge time	Time	90mins
Environment temperature	Degree C	+2-50
Atmospheric relative humidity		Less 80%
Adapter Import		110V-240V; 50-60Hz
Adapter output		3VDC, 0.5A
Dower dissipation		Less 2W

## Warranty specification

The company will undertake the exchange if the damaged which from material or manufacturing within one month from the date of delivery of this instrument, The company will provide the maintenance with maintenance of charge if the if the damaged which from material or manufacturing within 12 months from the date of delivery of this instrument, During the warranty period, the company will provide the free of charge repairing or replacement accessories if has been proved the instrument was damaged .

## Warranty coverage

If is misused 、 unsatisfactory operating and the damage caused by unauthorized repair or modification which shall not be covered by warranty .It appropriate the additional cost which repair outside the warranty period.

## Safety instructions

Please read this manual carefully and below safety instructions before using this instrument.

1. Only the original power adapter can be used for the instrument.
2. The instrument should be put in dry working environment.

## Packing List

Multi-channel motorize decapper	1 unit
USB cable	1 unit
Adapter	1 unit
Operation Manual	1 unit